


Dehydrogenation mechanism of LiBH₄ by Poly(methyl methacrylate)

Hydrogen is considered as a promising alternative energy carrier owing to its high-energy density, abundance, light weight and pollution-free burning [1]. Developing a safe and efficient hydrogen storage material is one of the key challenges for the mobile application of hydrogen [2]. Due to the high gravimetric (18.5 wt.%) and volumetric (121 kg H₂/m³) hydrogen density, lithium borohydride (LiBH₄) has been acknowledged as a potential candidate for hydrogen storage materials [3]. However, due to the unfavorable high thermal stability (e.g. decomposition peak temperature of ~ 470 °C), the practical utilization of LiBH₄ as hydrogen storage medium is hampered [4]. Hence, several approaches including reactant destabilization, catalyst/additive introduction, nanostructuring, and anion/cation substitution have been applied to decrease the dehydrogenation temperature and accelerate the kinetics [5].

Nanoengineering had been demonstrated to be a useful method to reduce the dehydrogenating/rehydrogenating temperature of LiBH₄ by decreasing diffusion path lengths and increasing surface areas [6]. However, nanoscale LiBH₄ is too reactive and very sensitive to the water and oxygen in the air, which impede its practical utilization. PMMA ((Poly (methyl methacrylate))), with a high permeability ratio of H₂/O₂, was reported to have good gas selectivity [7]. Therefore, in this project, PMMA was applied to protect LiBH₄ from oxygen and water but let the hydrogen get in or out freely (Scheme 1). Furthermore, the nanoconfinement of LiBH₄ in the fine network pore of PMMA and the interaction between the B atom in LiBH₄ and the O atom in C=O of PMMA resulted in a much lower hydrogen release temperature of LiBH₄. LiBH₄ PMMA composite started to dehydrogenate at 53°C and released 5.2 wt.% of hydrogen at 162°C within 1 h. This project provides a general strategy to utilize a gas-selective polymer to protect air-sensitive hydrogen storage compounds and improve their hydrogen storage properties.

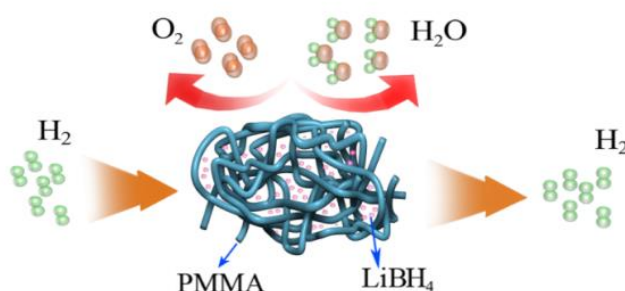

Scheme 1. Schematic illustration of LiBH₄ protected from oxygen and water by PMMA.

References

- [1] Jain, I. P. *Int J Hydrogen Energ* 2009, **34**, (17), 7368-7378.
- [2] Schlapbach, L.; Züttel, A. *Nature* 2001, **414**, (6861), 353-358.
- [3] Züttel, A.; Wenger, P.; Rentsch, S.; Sudan, P.; Mauron, P.; Emmenegger, C. *J Power Sources* 2003, **118**, (1-2), 1-7.
- [4] Saldan, I. *Cent Eur J Chem* 2011, **9**, (5), 761-775.
- [5] Li, H.-W.; Yan, Y.; Orimo, S.-i.; Züttel, A.; Jensen, C. M. *Energies* 2011, **4**, (1), 185-214.
- [6] Fang, Z. Z.; Wang, P.; Rufford, T. E.; Kang, X. D.; Lu, G. Q.; Cheng, H. M. *Acta Mater* 2008, **56**, (20), 6257-6263.
- [7] K. J. Jeon, H. R. Moon, A. M. Ruminski, B. Jiang, C. Kisielowski, R. Bardhan and J. J. Urban, *Nat. Mater.*, 2011, **10**, 286.

Project summary by:

Jianmei Huang, Liuzhang Ouyang, Min Zhu
School of Materials Science and Engineering
Key Laboratory of Advanced Energy Storage Materials of Guangdong Province,
South China University of Technology,
Guangzhou 510641
People's Republic of China

**Paper Reference:**

Jianmei Huang *et al.* (2015) "Dehydrogenation mechanism of LiBH_4 by Poly(methyl methacrylate)"
Journal of Alloys and Compounds **645** (1), S100-S102

Hidden Product:

[QGA](#)