

Absolute Quantification of Deuterium Ion Fluxes from an ECR Plasma Source

Plasma discharges are widely used in technical processes such as surface modification and thin-film deposition and etching. In this sense, the corresponding quantitative characterization of plasma sources is a topic of particular importance for a fundamental understanding of plasma sources as well as for a technological optimization of processes. Hydrogen plasmas play a certain role in technical discharges so that a precise determination of the absolute hydrogen fluxes to a plasma-treated surface is very desirable. This is important for understanding both the discharge mechanism and the plasma-induced processes at the surface.

A custom built electron-cyclotron-resonance (ECR) plasma source with a biased sample holder was thoroughly characterized with respect to the deuterium fluxes impinging on the substrate holder. The mass distribution of the impinging ions was measured with a HIDDEN plasma monitor (Model HIDDEN EQP 300). The plasma monitor was inserted in a mock-up sample holder thus allowing the measurement of the flux arriving at the usual sample position. In order to shield the plasma monitor from the magnetic field necessary for the ECR discharge, the mock-up sample holder was constructed from soft iron. An additional soft iron shielding cylinder was placed around the ion optics of the plasma monitor. The plasma monitor is differentially pumped in order to maintain a pressure of about 10^{-7} to 10^{-6} Pa during operation. The aperture in the mock-up sample holder was $10\mu\text{m}$. For absolute quantification of the mass-integrated ion flux to the sample holder a retarding field analyser (RFA) was used.

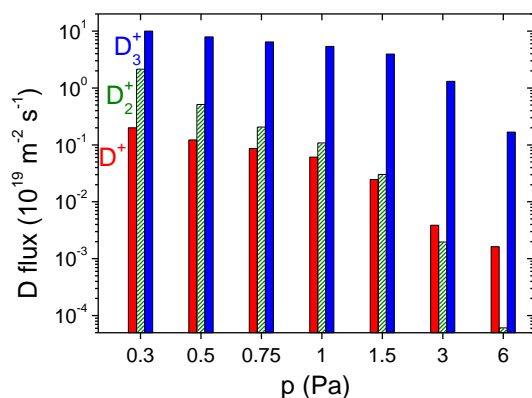


Figure 1: Ion-species-resolved deuterium flux as a function of the D_2 gas pressure, measured at a constant microwave input power of 144 W and with the sample holder at floating potential

The resulting ion fluxes impinging on the substrate holder for a microwave power of 144 W are presented in Fig. 1 as a function of operating pressure in the range from 0.3 to 6 Pa. Even at the lowest investigated pressure of 0.3 Pa the dominating ion species hitting the sample surface is D_3^+ . The dominant production pathway for this ion is an ion-molecule reaction in the bulk plasma. A D_2^+ ion collides with a D_2 molecule producing a D_3^+ and a D atom. With increasing pressure the probability for this reaction increases so that the relative ion species composition changes. With increasing pressure more and more D_3^+ ions are produced at the expense of the primary D_2^+ ion which is produced in the bulk plasma by electron-induced ionization of D_2 molecules. At pressures higher than about 1.5 Pa the total target current decreases due to a decrease of the plasma density, but still the relative

contribution of D_3^+ continues to increase. At the highest investigated pressure of 6 Pa the contribution of D_2^+ is almost negligible (less than 10^{-3}). In contrast to D_2^+ the fraction of D^+ does not significantly change. D^+ is, as D_2^+ , produced by electron-induced ionization of D_2 molecules and additionally by direct ionisation of atomic D (D plasma concentration is about 1 %), but in contrast, the cross section for an ion molecule reaction of D^+ is much lower than that of D_2^+ .

For the defined standard operating parameters of our plasma source, i.e. a D_2 plasma at $p = 1.0$ Pa, $P_{MW} = 144$ W and the sample holder at floating potential, the total deuteron flux in the form of ions is $5.6 \times 10^{19} \text{ Dm}^{-2}$. This corresponds to a total ion flux of $1.9 \times 10^{19} \text{ ions m}^{-2}$. 94% of these ions are D^{+3} , 3% are D^{+2} , and 3% are D^+ ions. The ion flux increases monotonically by about a factor of 2 if the DC bias voltage is increased to -600 V.

Experimental setup:

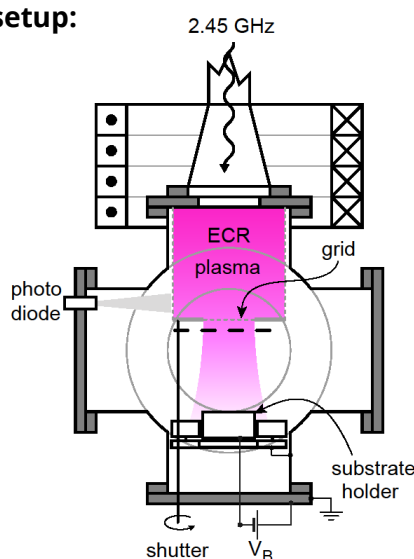


Figure 2: Schematic view of the experimental set-up. To quantify the particle flux to the substrate the substrate holder is replaced by a mock-up containing the EQP plasma monitor (not shown)

Project Summary by:

A. Manhard, T. Schwarz-Selinger, and W. Jacob
 Max-Planck-Institute für Plasmaphysik, EURATOM Association,
 Boltzmannstr. 2, 85748 Garching, Germany



Reference:

A. Manhard, T. Schwarz-Selinger, and W. Jacob (2011) "Quantification of the Deuterium Ion Fluxes from a Plasma Source" *Plasma Sources Science and Technology* **20** 015010
 doi:10.1088/0963-0252/20/1/015010

Paper Reference:

P. Wang, W. Jacob, M. Balden, A. Manhard, T. Schwarz-Selinger (2012) "Erosion of tungsten-doped amorphous carbon films in oxygen plasma" *Journal of Nuclear Materials* **420** (1-3), 101-109

Hidden Product:

EQP Mass and Energy Analyser

Follow the link to the product catalogue on our website for further information

<http://www.hiddenanalytical.com/en/products/for-thin-films-plasma-surface-engineering/eqp>