


Hiden HPR-60


Molecular Beam Mass Spectrometer (MBMS) for the quantitative analysis of reactive gas species

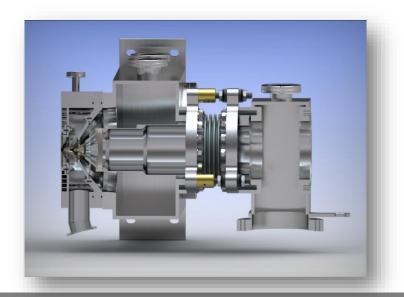




HPR-60 Overview

- The Hiden HPR60 Molecular Beam Mass Spectrometer (MBMS) is a compact, mobile gas analysis system for the quantitative analysis of reactive gas species.
- Radicals, ions, polymers and clusters are sampled via a multistage differentially pumped inlet, forming a molecular beam that is transferred to the ion source of a precision triple filter quadrupole mass spectrometer.

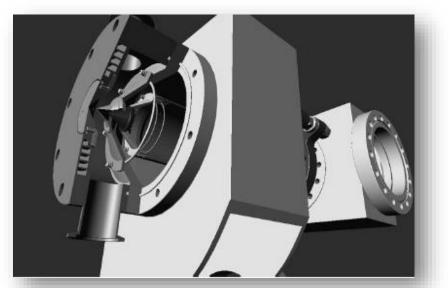


HPR-60 System Schematic

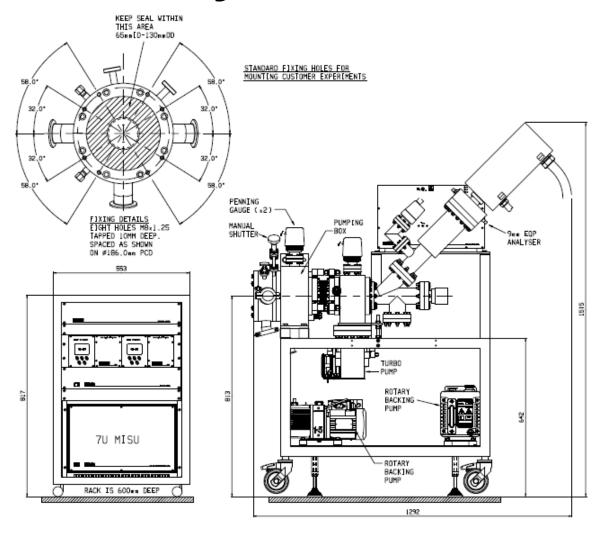

HPR-60 Features

- Molecular beam sampling at atmospheric pressure
- +ve and -ve ion analysis
- User replaceable skimmer
- Cones (can be biased)
- Electron attachment ionisation mode for the study of electro-negative radicals
- APSI-MS soft ionisation mode for radicals analysis
- Mass range options: 300, 510 or 1000 amu
- Energy range options 100 eV or 1000 eV



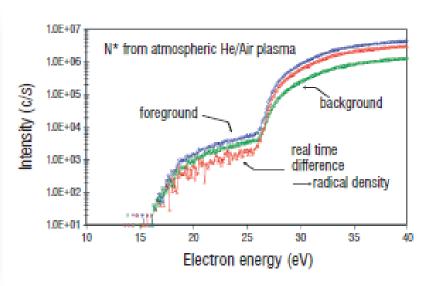


HPR-60 Applications

Both two and three stage differentially pumped versions are available to address a broad range of applications covering the pressure range 10-4 mbar to atmospheric, including reaction kinetics in;

- Environmental and atmospheric chemistry
- Low and high pressure plasma chemistry
- Catalytic reactors
- CVD / MOCVD
- Combustion chemistry
- Flame chemistry




HPR60 Typical Operating Pressures

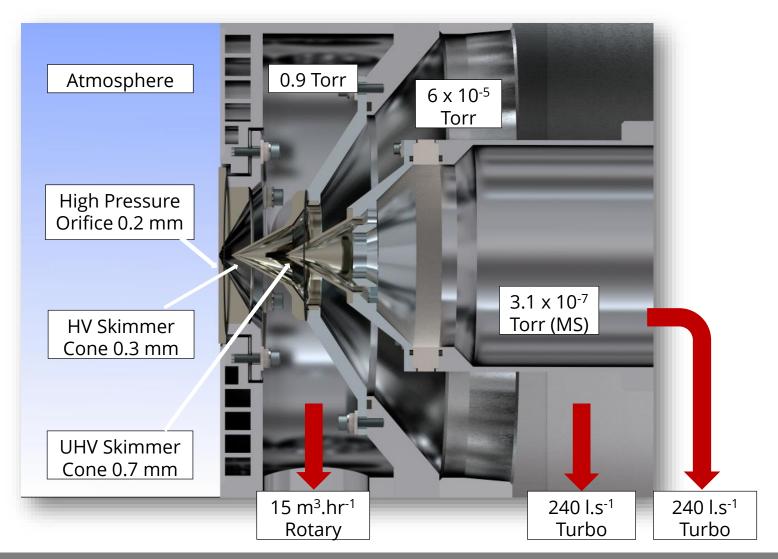

HPR-60 Water Cooled Stage for High Temperature Measurements


- Water cooled channels are offered as standard, with optional chiller.
- Special graphite seal allows heat conduction from the orifice.
- Assuming water cooling is used at a flow rate of ~1 l.min⁻¹, with 20 °K rise. A cooling power of 1.3 kW is obtained.


For higher temperature applications, ceramic cones and heat shield are also offered.

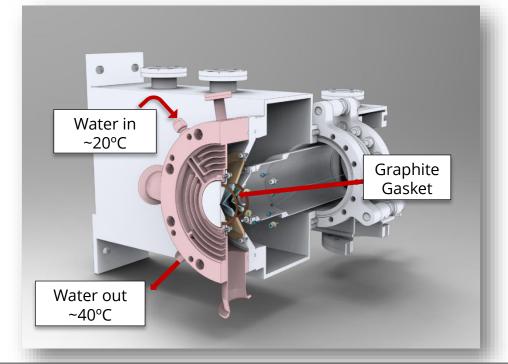

HPR-60 Cones

- Control the pressure differential obtained through stages 1, 2, and 3.
- Standard orifice sizes– suitable for sampling from around atmospheric pressure as standard. Other pressure regimes are catered for.
- Radicals are sampled via a multi stage, differentially pumped skimmer inlet. Sampling is carried out in the supersonic 'mach disc' region, interactions with other species and chamber walls are negligible.
- Other metal and ceramic cones available Please contact Hiden for more details.

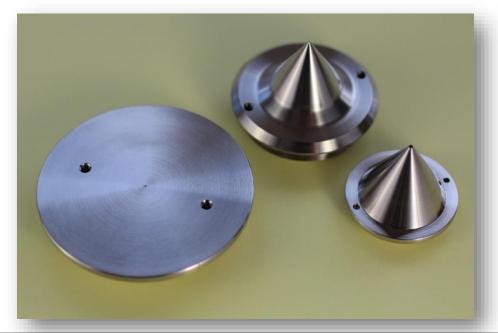

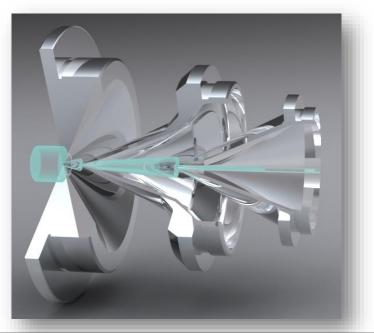

HPR-60 Mass Spectrometer Selection

EPIC Probe

- Sub PPM detection of plasma ions, neutrals and radicals.
- Mass ranges of 0-50, 0-300, 0-500
 & 0-1000 amu available.

EQP Probe


- Sub PPM detection of plasma ions, neutrals and radicals.
- Ion Energy Analysis, 45° sector energy analyser with 0.25 eV resolution
- Mass ranges of 0-50, 0-300, 0-500
 & 0-1000 amu available.



EPIC Probe – Mass Analyser



EQP Probe – Mass/Energy Analyser

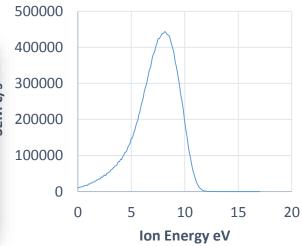


Typical EQP Energy Spectrum from an atmospheric Plasma Discharge.

EQP Extraction Optics



- Software controlled Extraction / Focus
- Discriminates +ve, -ve, e-, radicals
- Fully tuneable for optimal detection
- Integrated ionisation source

EQP 45° Electrostatic Sector Analyser

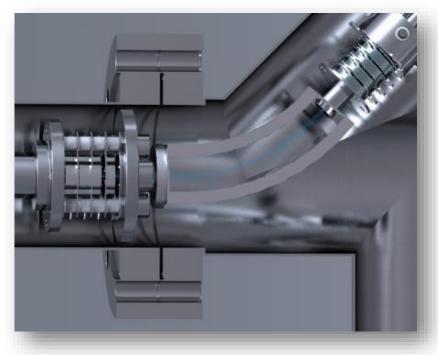

- Optimum transmission cf. Bessel box.
- Minimum perturbation of ion flight path.
- Constant ion transmission at all energies.
- Energy resolution down to 0.25 eV FWHM.
- Energy scan at increments from 0.05 eV .
- Floating option up to 10 KeV.


Triple Filter Mass Spectrometer

Why have a triple filter?

Two main advantages:

- Strict control over the quadrupole entrance and exit fields provides enhanced sensitivity for high mass transmission and increased abundance sensitivity
- 2. Enhanced long-term stability. The bulk of the deselected ions from the quadrupole ioniser deposit harmlessly on the RF-only pre-filter stage, minimising contamination on the mass selective primary filter


Secondary Electron Multiplier (SEM) Detector

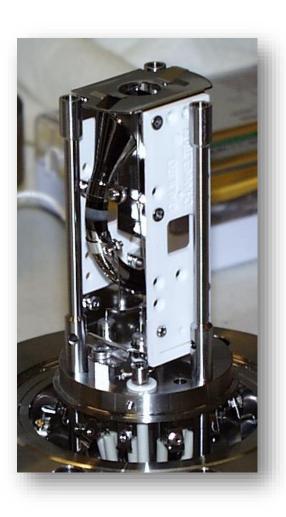

- 7 decade continuous dynamic range.
- 24 bit counter for 1c /s resolution.
- Faraday Cup option for high density plasmas.
- Signal gating with 1 μs resolution energy & mass distributions vs time.
- TTL / DDE data export options.

Summary

- Molecular beam Mass Spectrometer (MBMS)
- Designed and manufactured by Hiden in the UK.
- Radicals, ions, polymers and clusters are sampled via a multistage differentially pumped inlet, forming a molecular beam that is transferred to the ion source of a precision triple filter Quadrupole mass spectrometer.

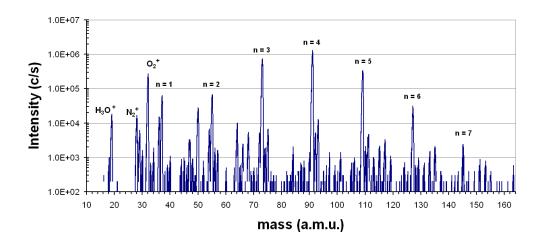



Selected Publications

- Ambient air particle transport into the effluent of a cold atmosphericpressure argon plasma jet investigated by molecular beam mass spectrometry. 2013. M Dünnbier et al. *J. Phys. D: Appl. Phys.* **46** 435203
- The reaction mechanism of the spray Ion Layer Gas Reaction process to deposit In₂S₃ thin films. 2011. S Gledhill et al. *Thin Solid Films* 519 6413-6419
- Atmospheric pressure plasma analysis by modulated molecular beam mass spectrometry. 2006. Y Aranda Gonzalvo et al. *J. Vac. Sci. Technol.* A 24(3) May/June
- A plasma needle generates nitric oxide. 2006. E stoffels et al. *Plasma Sources Sci. Technol.* **15** 501-506
- A mass spectrometric study of ions extracted from a point-to-plane dc corona discharge in N₂O at atmospheric pressure. 2008. JD Skalny et al. J.Phys. D: Appl. Phys. 41 085202
- Positive Ion Mass Spectrometry during an Atmospheric Pressure Plasma Treatment of Polymers. 2009. AJ Beck et al. *Plasma Process.* Polym. 6 521-529







Hiden HPR-60 Users

- Samsung Electronics
- Xian Jiaotong University
 - KAUST
- Old Dominion University
- University College London
- Institute of Plasma Physics
- Atmospheric Afterglow Technologies
 - University of Liverpool
 - CEA LETI
 - CBTE
 - Nagoya Institute
 - Oak Ridge National Laboratory
 - Paul Scherrer Institut











Quadrupole Mass Spectrometers for Advanced Science



Hiden Analytical Ltd. 420 Europa Boulevard Warrington, WA5 7UN, England

www.HidenAnalytical.com

info@hiden.co.uk

Tel: +44 (0)1925 445 225

