

Download procedure

The following instructions apply to Microsoft Edge, other web browsers can be used and have equivalent dialog boxes.

1. Enter the address
https://ftp.hiden.co.uk
into your web browser's address bar.
2. Press the **Enter** key.

If you are viewing this document in an electronic format you may be able to copy and paste the address. Usually, using **Ctrl+C** and **Ctrl+V** will respectively copy and paste.

The **Login** page, as shown in Figure 1, will be displayed.

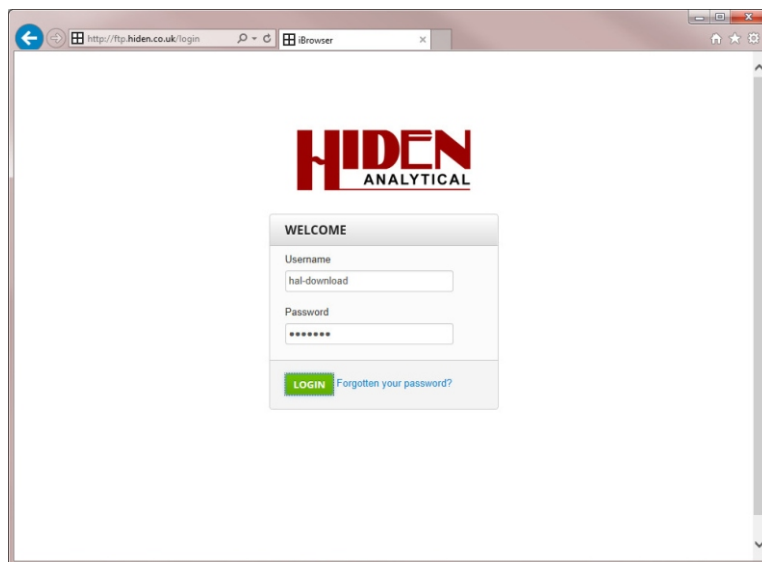


Figure 1 The Login page

3. In the box labelled **Username**
type **hal-download**
4. In the box labelled **Password**
type: **hiden07**
5. Click the **Login** button.

The Username and Password are case sensitive.

The web browser will display the **Spaces** page, see Figure 2.

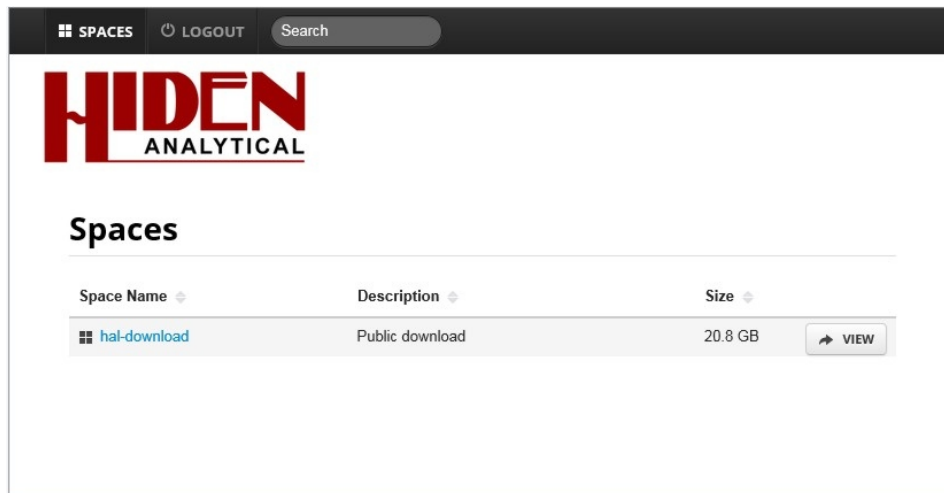


Figure 2 The Spaces page

6. Click the **hal-download** link.

The **hal-download** page will be displayed. See Figure 3.

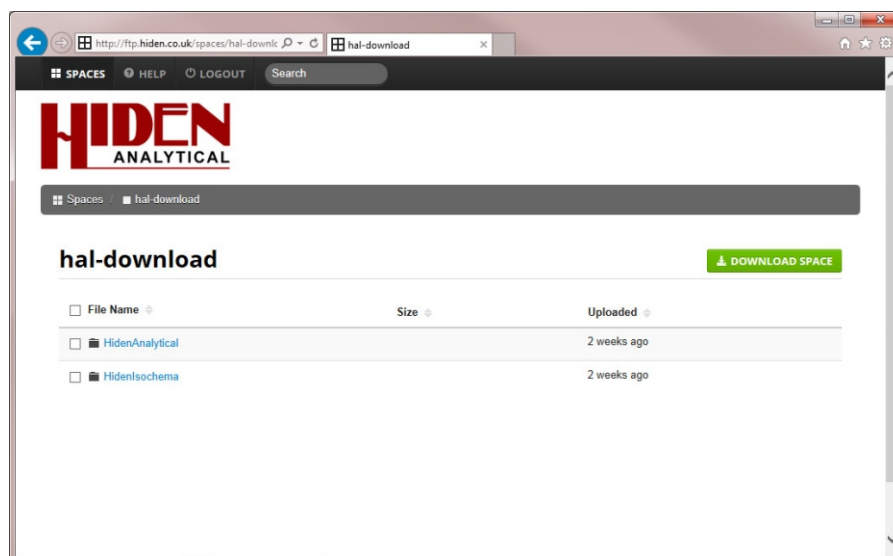


Figure 3 The hal-download page

7. Click the **HidenAnalytical** link.

The HidenAnalytical options will be displayed. See Figure 4.

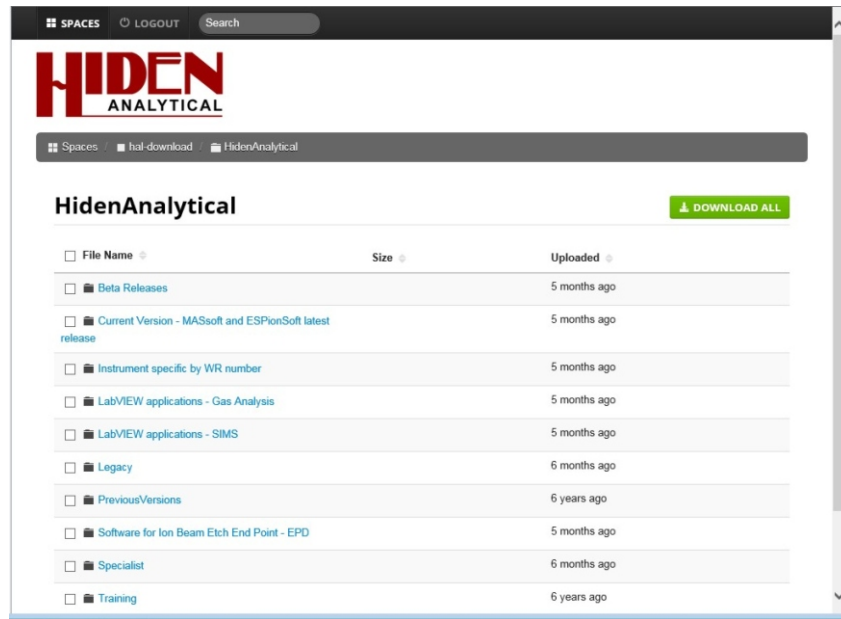


Figure 4 The HidenAnalytical options page

Note

The exact directory structure of the FTP site may change from time to time.

8. Click on the **CurrentVersion-MASsoft and ESPionSoft latest release** link.

The Current Version page is displayed as shown in Figure 5.

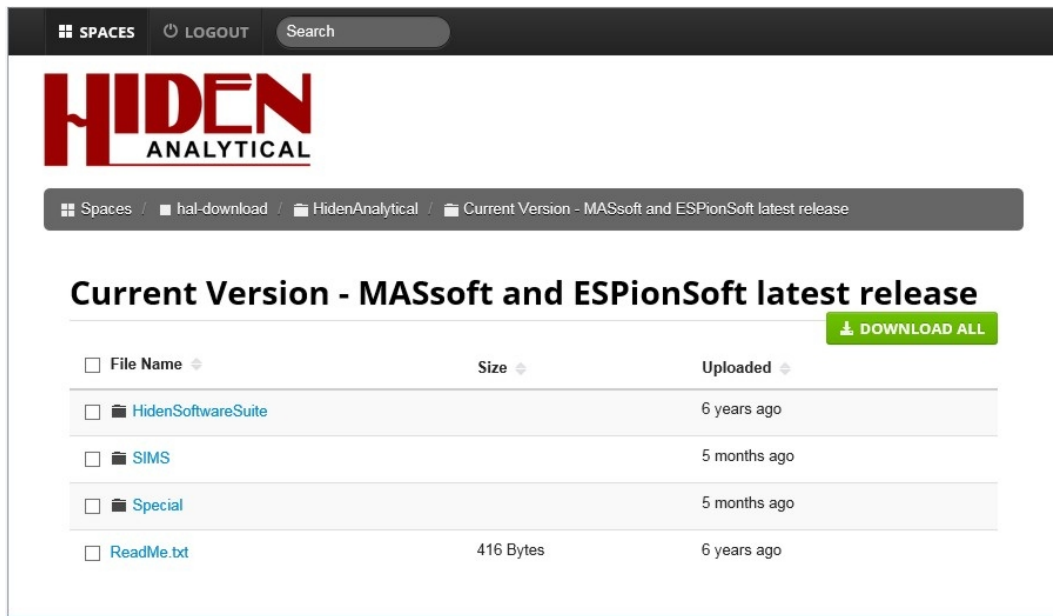


Figure 5 Current Verison page

9. Click on the **HidenSoftwareSuite** link.

The HidenSoftwareSuite page is displayed as shown in Figure 6.

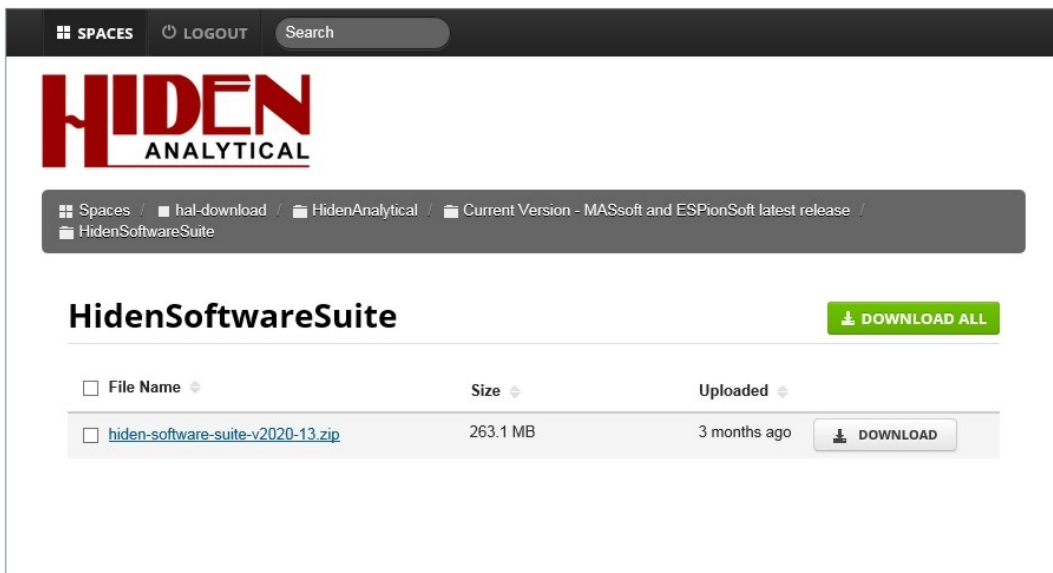


Figure 6 HidenSoftwareSuite page

- | | |
|--|--|
| 10. Click on the DOWNLOAD button next to the hiden-software-suite-v202x-xx.zip link. | The file will begin downloading automatically. |
| 11. When the download is complete right-click on the file and select Show in folder . | Windows Explorer will be displayed. |
| 12. Right-click on the hiden-software-suite-v202x-xx.zip file. | A pop-up menu will be displayed. |
| 13. In the pop-up menu click on Extract All... | The Extract Compressed (Zipped) Folders dialog box will be displayed, as shown in Figure 7. |

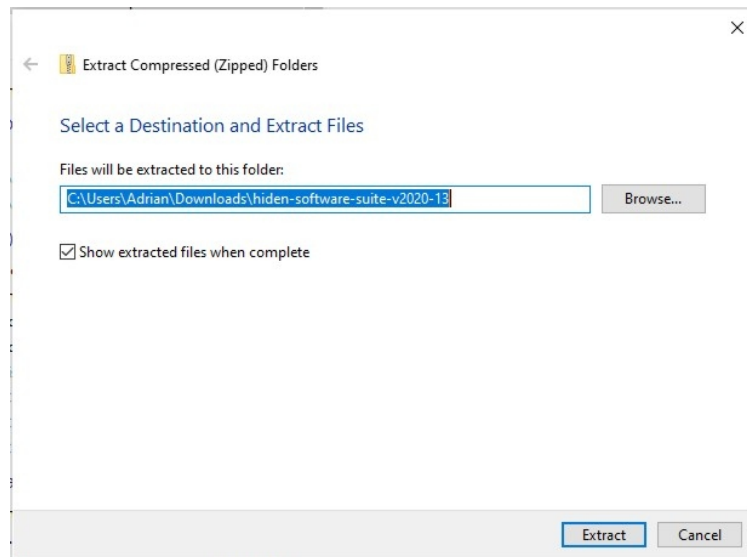


Figure 7 Extract Folders dialog box

- | | |
|--|--|
| 14. Click on the Browse... button. | The Select a destination dialog box will be displayed. |
| 15. Select a folder to extract the files to then click OK . | The dialog box will close. |
| 16. Click Extract . | The Copying dialog box will be displayed, as shown in Figure 8. |

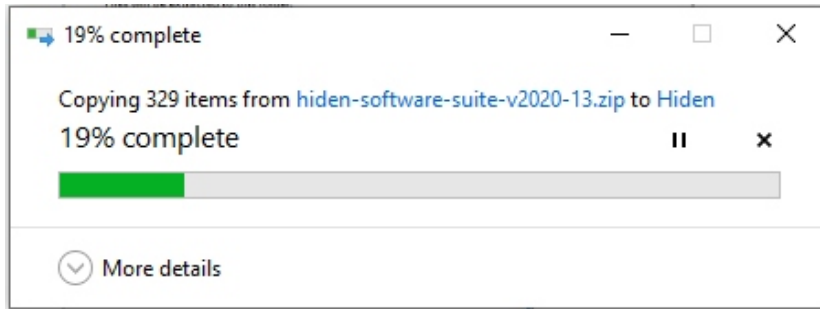


Figure 8 Copying files dialog box

17. When copying is complete navigate to the folder specified in step 15.
18. Double click on the folder named **hidden-software-suite-v202x.xx.zip**. A list of files similar to the one shown in Figure 9 will be displayed.

Note

The **hidden-software-suite-v202x-xx** folder contains a file called **issue.txt** which lists all of the improvements made to the software since the last release.

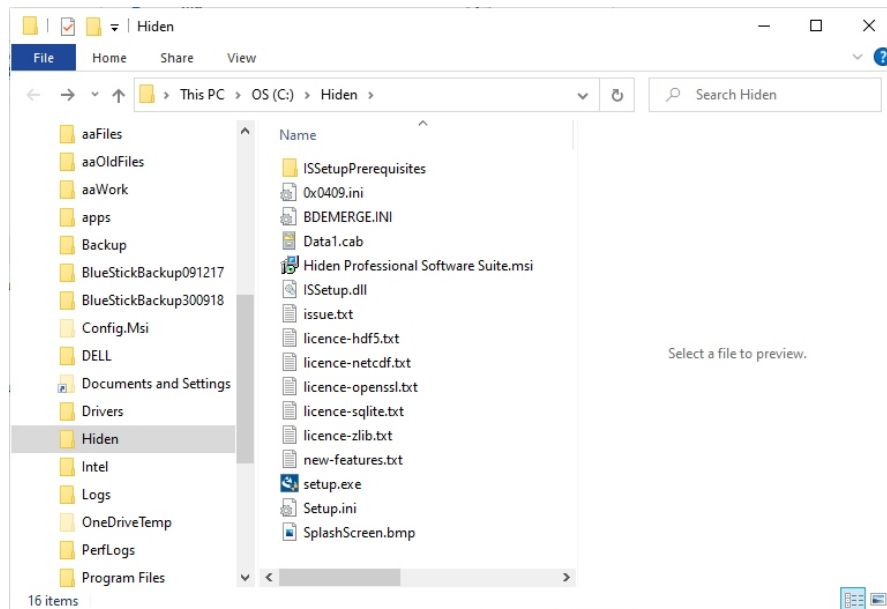


Figure 9 hidden-software-suite-v202x-xx in Windows Explorer

19. Double-click on the **setup.exe** file.

The **User Account Control** dialog will be displayed.

20. Click the **Yes** button.

The software installation will begin.