

#### **Hiden HPR-20 TMS**

Transient MS for Fast Event Gas Analysis







#### Introduction

The Hiden HPR-20 TMS is configured for continuous analysis of gases and vapours at pressures near atmosphere.

Operating to 200°C, the QIC (quartz inert capillary) flexible 1m capillary inlet provides fast response times of less than 150 ms.

The HPR-20 TMS system has a mass range of 200 amu (300, 510, 1000 amu options) and a detection capability from 100% to less than 5 ppb.

The TMS system is optimised for fast response studies and can respond to changes in gas composition with a 5 decade response time in <200 ms.



#### TMS – for Fast Event Studies

Features of the system to optimise the response for fast event studies include:

- Pulse Ion Counting detection with 7 decade continuous log scale
- 1m capillary gives <150 response time
- Open ion source increases pumping speed on ion source





# **Ultra-fast response**



HPR-20 TMS: > 5 decades response in < 0.2 s




# **Ultra-fast response**



The HPR-20 TMS tracks the changes in concentration of fast gas pulses with incredible response and accuracy over more than 5 decades. This data shows the measurement of 5 pulses in 5 seconds.



#### **HPR-20 TMS Vacuum Schematic**

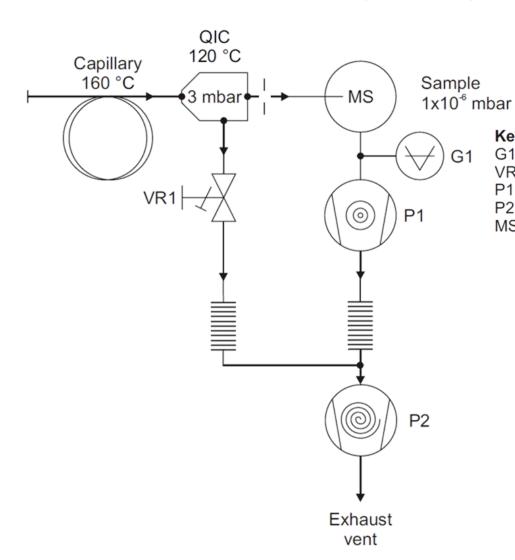

#### Key

G1 Penning gauge VR1 QIC Inlet bypass control valve P1 60 l/s turbo drag pump P2 Backing and bypass Scroll pump MS UHV Housing (Mass spectrometer chamber)




Backing and bypass Scroll Pump



# **HPR-20 TMS Analyser:**

#### **Hiden HAL 3F PIC Triple Filter Mass Spectrometer**

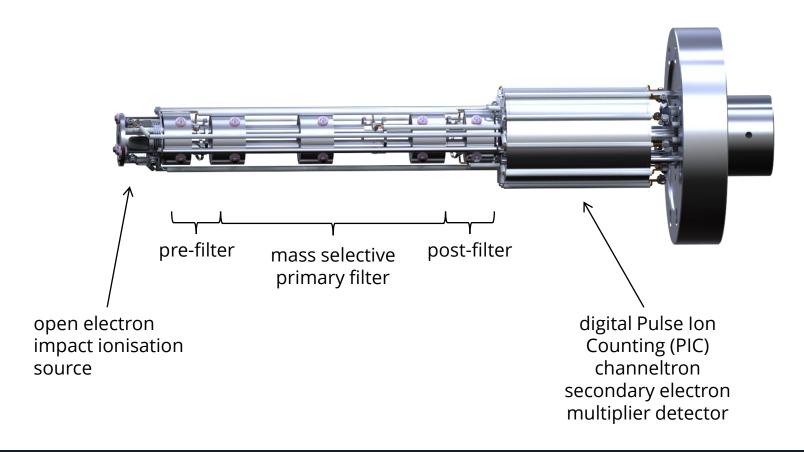



# **Triple Filter Mass Spectrometer**


Why have a triple filter?

Two main advantages:

- 1. Strict control over the quadrupole entrance and exit fields provides enhanced sensitivity for high mass transmission and increased abundance sensitivity
- 2. Enhanced long-term stability. The bulk of the deselected ions from the quadrupole ioniser deposit harmlessly on the RF-only pre-filter stage, minimising contamination on the mass selective primary filter.



# **QIC Inlet Technology**

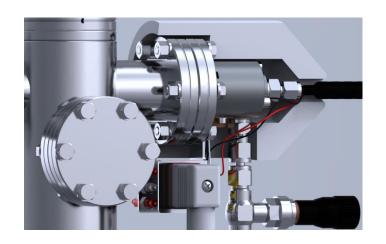

Quartz and Platinum Wetted Surfaces No memory effects

Heated Capillary No condensation effects


Flow Matched → Optimum response / recovery

Minimal Internal Volume PPB detection

Interchangeable Sampling Capillaries —— Analysis from 10 mbar to 2 Bar



# **Typical Mass Spectrum of Air**

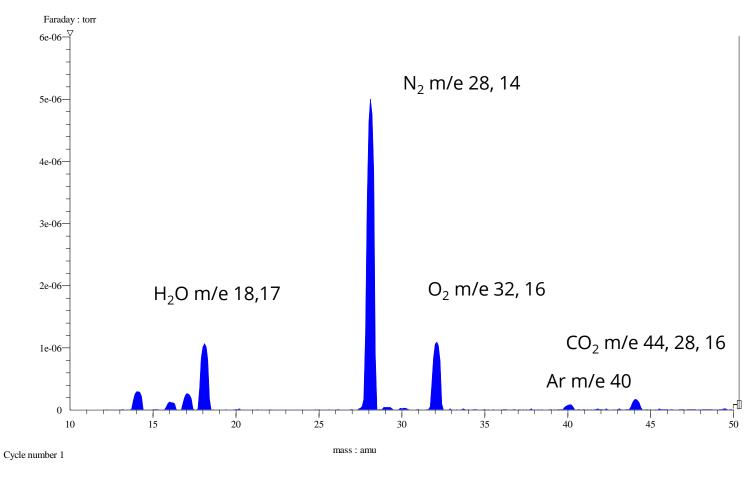

Note: Different species can have the same mass e.g. CO,  $N_2$  m/e 28



#### **Soft Ionisation**

Unique to Hiden gas analysis systems, soft ionisation allows users to selectively ionise different gases by setting the ionisation energy for a particular mass.

This powerful technique can simplify the analysis of otherwise complex cracking patterns from multi-component gas/vapour mixtures.

The ionisation energy can be altered from 4 to 150 eV, in 0.1 eV increments. Standard operation is at 70 eV.


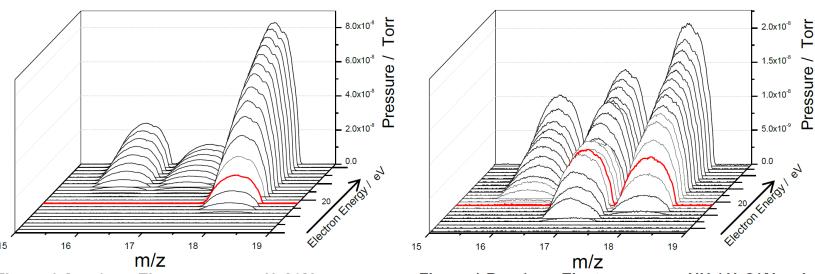


Figure 1 A: m/z vs Electron energy-H<sub>2</sub>O/Air

Figure 1 B: m/z vs Electron energy-NH<sub>3</sub>/ H<sub>2</sub>O/Air mix



#### **MASsoft Professional control software**

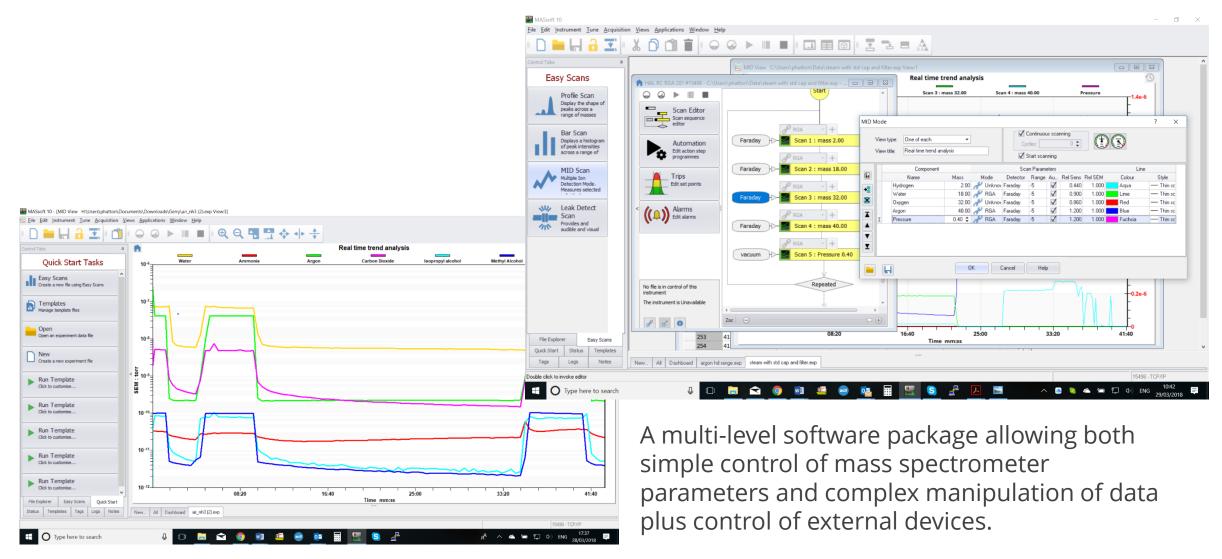



## **HPR-20 TMS - Applications**

- Fast event studies
- Reaction kinetics
- Selective Catalytic Reduction (SCR)
- Steady State Isotopic Transient Kinetic Analysis (SSITKA)
- High speed switching analysis
- Operando studies
- SpaciMS





#### **Applications: Catalysis Research**

A kinetic study of the effect of  $H_2$  on the Selective Catalytic Reduction of  $NO_x$  with octane using isotopically labelled <sup>15</sup>NO, using an HPR-20 QIC TMS system.

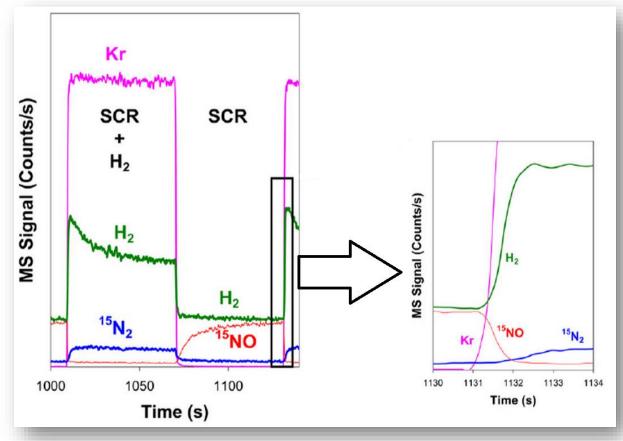



Figure 1. Mass spectrometer response for <sup>15</sup>N<sub>2</sub>, <sup>15</sup>NO, H<sub>2</sub> and Kr when switching 0.72% H<sub>2</sub> in and out of a SCR feed stream over the catalyst at 300°C.

Ref: J. P. Breen, R. Burch, C. Hardacre, C. J. Hill and C. Rioche (2007). J. Catal., 2007, 246, 1, p1-9.



# **Applications: Breath Analysis**

Breath by Breath Analysis of Expired Isoprene during Exercise

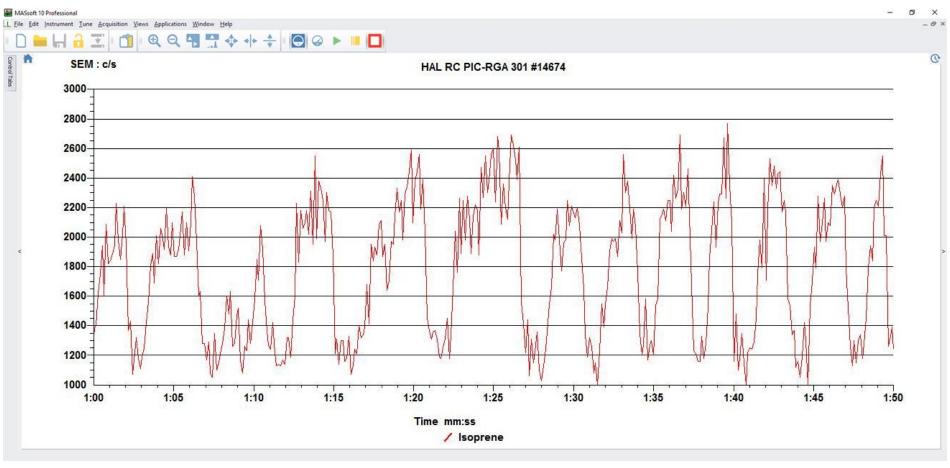

HPR-20 TMS data showing breath by breath isoprene levels during an exercise test.












#### **Imperial College** London

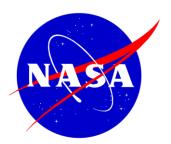





Eidgenössische Technische Hochschule Zürich Swiss Federal Institute of Technology Zurich

#### **Hiden HPR-20 Users**

NASA Dow Chemical Exxon-Mobil Imperial College MIT

University of British Columbia University of Queensland **BASF** 

Seoul National University Suzuki

University of Cambridge Beijing Institute of Technology

Samsung ETH Zürich **KAUST** 

**Durham University** Siemens Shell







**Massachusetts Institute of Technology** 















SAMSUNG





## **Summary**

- Specifically designed for fast event studies less than 150 ms response time
- Bench-top triple filter quadrupole mass spectrometer gas analysis system
- Real-time, multi-species analysis 5 PPB to 100%
- Soft ionisation for reduced spectral fragmentation and simplified data interpretation





