
www.hidenanalytical.com info@hiden.co.uk

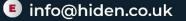

Hiden QGA

Quantitative Gas Analysis System







Hiden QGA System



A compact bench-top gas analyser configured for real time analysis.

www.HidenAnalytical.com 📧 info@hiden.co.uk



Introduction


The Hiden QGA quantitative gas analysis system is configured for continuous analysis of gases and vapours at pressures near atmosphere.

Operating to 200°C, the QIC (quartz inert capillary) flexible 2m capillary inlet provides fast response times of less than 300ms.

The QGA system has a mass range of 200 amu (300 amu option) and a detection capability from 100% to less than 100ppb.



QGA Vacuum Schematic

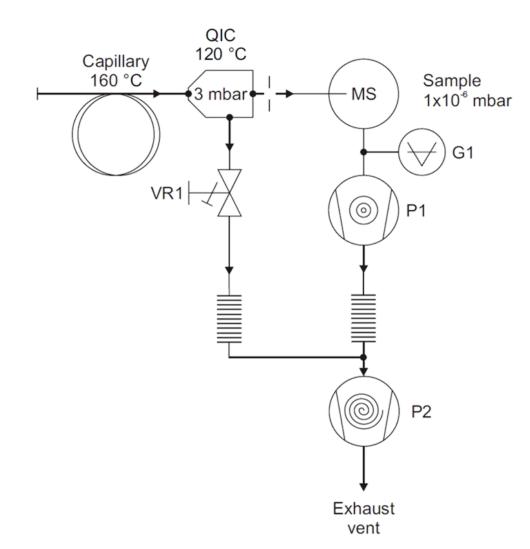

Key

토 info@hiden.co.uk


G1 Penning gauge VR1 QIC Inlet bypass control valve P1 60 I/s turbo drag pump P2 Backing and bypass Scroll pump MS UHV Housing (Mass spectrometer chamber)



Backing and bypass Scroll Pump




QGA Mass Spectrometer

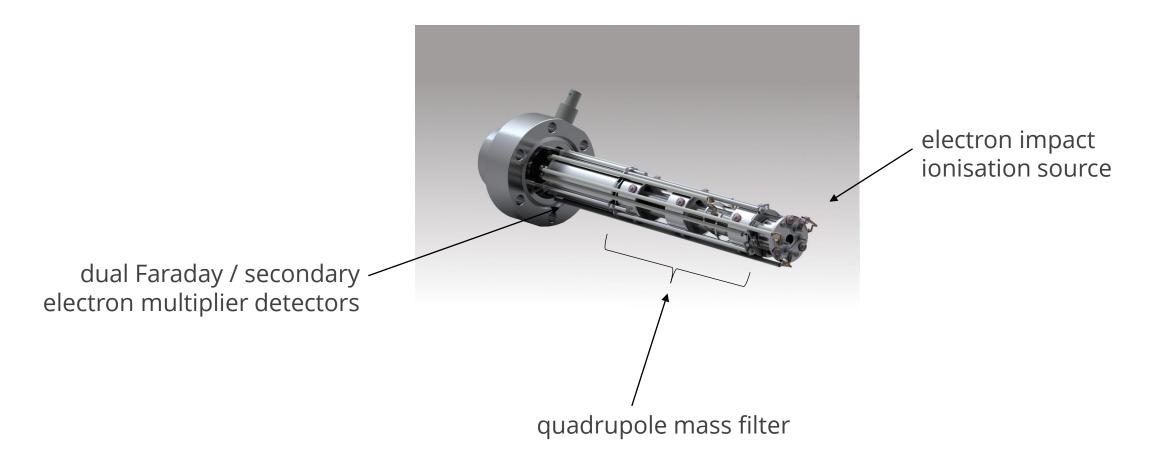




QIC Inlet Technology

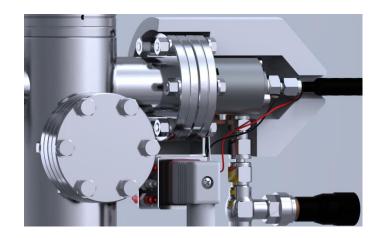

- Quartz and Platinum Wetted SurfacesNo memory effectsHeated CapillaryNo condensation effectsFlow MatchedOptimum response / recoveryMinimal Internal VolumePPB detectionInterchangeable Sampling CapillariesAnalysis from 10 Torr to 2 Bar
 - www.HidenAnalytical.com 💿 info@hiden.co.uk




Fast Response to Permanent Gases / Vapours

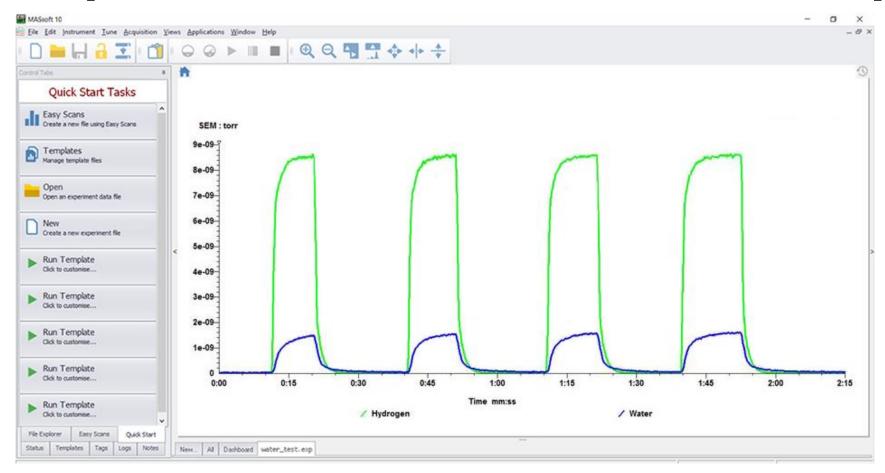

Data shows the response of a QGA system to gas and vapour during switching between a dry He stream and a wet H_2 and Ar flow. For clarity, only the H_2 and H_2O data is shown in the graph.



Typical Mass Spectrum of Air

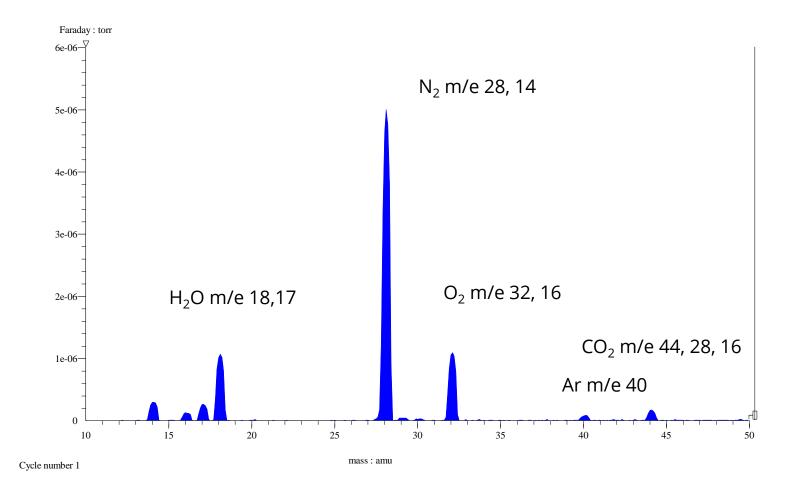

Note: Different species can have the same mass e.g. CO, N_2 m/e 28

www.HidenAnalytical.com 💿 info@hiden.co.uk




Soft Ionisation

Unique to Hiden gas analysis systems, soft ionisation allows users to selectively ionise different gases by setting the ionisation energy for a particular mass.

This powerful technique can simplify the analysis of otherwise complex cracking patterns from multi-component gas/vapour mixtures.

The ionisation energy can be altered from 4 to 150 eV, in 0.1 eV increments. Standard operation is at 70 e'

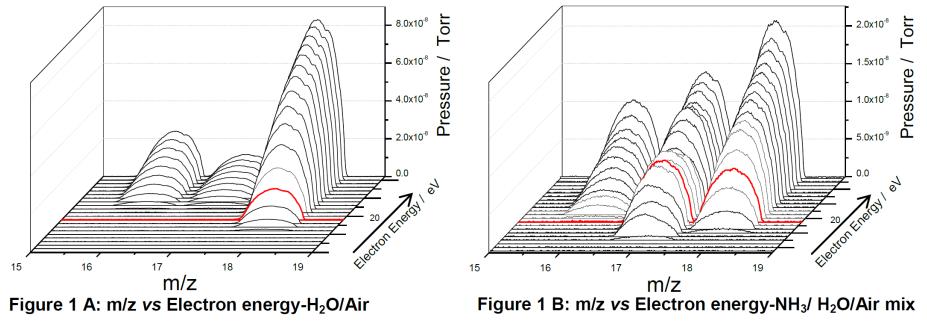




Gas Sampling Options

Pressure:

Inlet options are available for sampling both above and below atmospheric pressure. High pressure inlets for sampling at up to **30** bar and special capillaries for sampling down to **1** mbar.

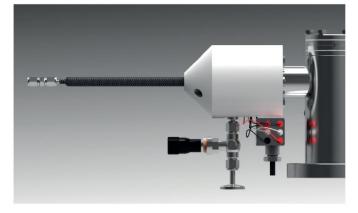





Multi-stream selectors:

2, 8, 16, 20, 40 and 80 way options

Temperature:

- Heated capillary extensions
- High temperature capillary inlets
- Hot-zone adaptors
- Heated multi-stream inlets



Type here to search

📄 숙 🧔


1 🗆

w] 🖆 📾

o 🗹

S

MASsoft Professional Control software

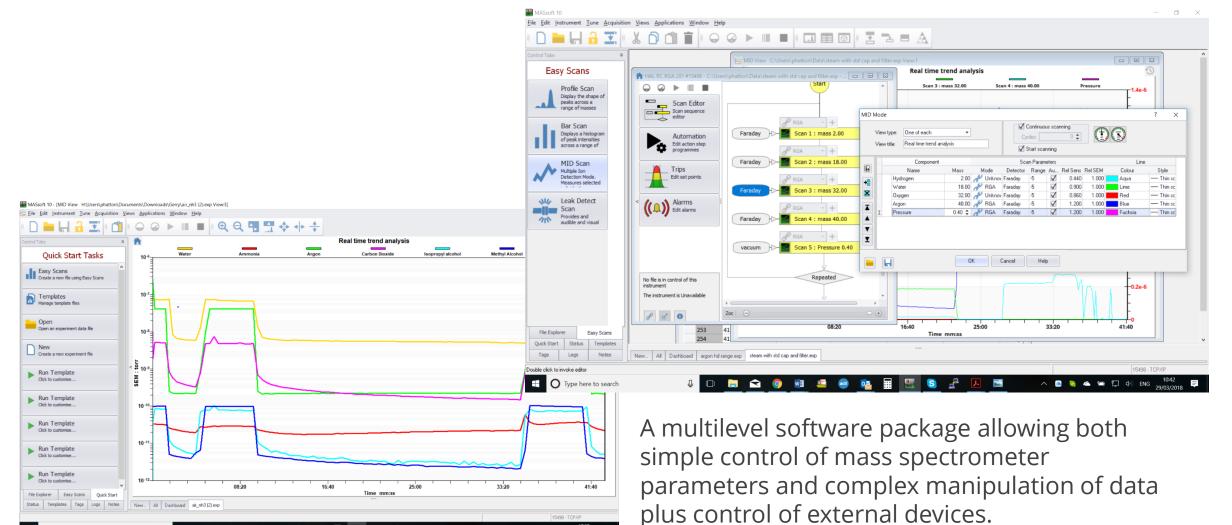

📌 🔨 📥 🚍 🖽 ENG



QGA Professional Software for Quantitative Gas Analysis



An application specific software package for quantitative gas and vapour analysis providing real time continuous analysis of up to 32 species with concentrations measured in the range 0.1PPM to 100%.


- Automatic subtraction of spectral overlaps
- Automated calibration routines
- Mass spectral library with intelligent scan feature
- Multi-stream support

www.hidenanalytical.com info@hiden.co.uk



QGA Professional Software for Quantitative gas Analysis



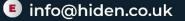



Applications


- Catalysis
- Reaction Kinetics
- TPD/TPR/TPO
- Thermal Analysis Mass Spectrometry
- Gas Purity Analysis
- Process Characterisation
- Fermentation Off Gas Analysis
- Environmental Gas Analysis
- Combustion Studies
- CVD/MOCVD





Applications: Catalysis Research

- Catalyst characterisation
- Kinetic and thermodynamic measurements
- TPD, TPO, TPR, TP-Reaction
- On-line continuous product analysis
- Total Surface Area / Metal Surface Area
- Mechanisms of Surface Reactions
- Heats of Adsorption and Coadsorption
- Operando Studies

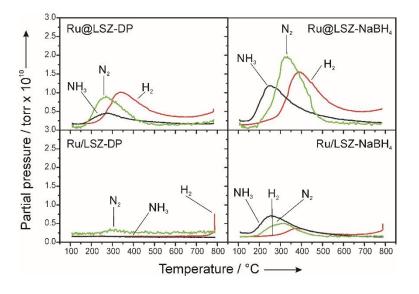


NH₃-TPD on embedded Ru@LSZ and impregnated Ru/LSZ catalysts.

Ref: Lorenzut et al. (2011) *Hydrogen production through alcohol steam reforming on Cu/ZnO-based catalysts*, Applied Catalysis B, **101** (3&4), 397-408.




Applications: Environmental Gas Analysis



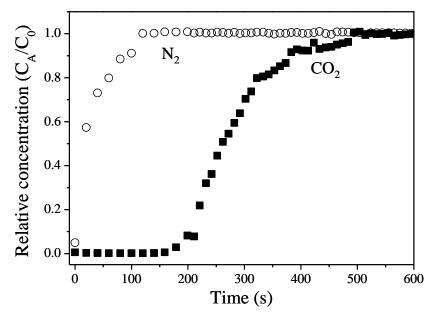

Breakthrough curve of CO_2 (15% CO_2 , 85% balance N_2) on mesoporous alumina. Ref: Yang et al. (2010) CO_2 adsorption over ionexchanged zeolite beta with alkali and alkaline earth metal ions, Mesoporous Materials **135** (1-3), 90-94.



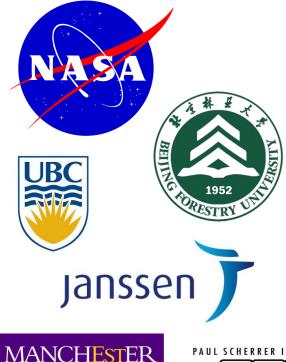

PAUL SCHERRER INSTITUT

The University of Manchester

1824

Imperial College London



Delft University of Technology

Hiden QGA Users

- Johnson Matthey
- Imperial College London
 - NASA
 - Poitiers University •
 - Nissan
 - **KTH Stockholm**
- Karlsruhe Institute of Technology
- University of Florida
- The University of Hong Kong
- Paul Scherrer Institut
- ULB Brussels University
- Texas A&M University
- University of Sao Paolo
 - University College London
- Technical University Denmark

- **Beijing Forestry University**
- CSIR Indian Institute of Petroleum
- Diamond Light Source ٠
- Shanghai Institute of Technology
- Janssen Pharmaceutica
 - PDVSA
 - Air Liquide
 - BASF
 - TU Delft
- Seoul National University
- University of Manchester
 - University of British Columbia
 - USTB
- University of Nottingham















Summary

- Compact bench-top gas analysis system
- Real-time, multi-species analysis 100 ppb to 100%
- Fast response to permanent gases and vapours less than 300 ms response time
- Automated calibration routines for accurate quantitative analysis
- Soft ionisation for reduced spectral fragmentation and simplified data interpretation